



## **INSTRUCTION FOR MIN MAX THERMOMETER**

- Contents:**
- 1 Thermometer**
  - 1 Battery**
  - 1 Cord**
- 
- Step 1**                    **Insert battery.**
- Step 2**                    **Attach cord to Min Max.**
- Step 3**                    **Feed cord into fridge via the hinged part of the door opening.**
- Step 4**                    **Insert the probe into a vaccine box and tape securely to middle of middle shelf.**
- Step 5**                    **Tape the cord to the shelf and to the side of the fridge being careful not to locate it near the hinged area of the door (this may cut the cord). Secure cord on the outside of the fridge and display the Min Max on top of the fridge for reading purposes.**
- Step 6**                    **The Min Max may be moved from top to bottom shelves each 24 hours to assess the temperature variations within the fridge.**
- Step 7**                    **Readings are done 2 x daily – am/pm.**
- Step 8**                    **Reset the Min Max after each reading by pressing the reset button.**
- Step 9**                    **The Min Max displays the current temperatures and also the minimum and maximum temperatures.**
- Step 10**                    **Please remove battery from Min Max and replace in plastic bag prior to returning to Tasmania Medicare Local South.**